

Decipher Mechanisms of Gene Expression Changes with Single-Cell Multiomics

ResolveOME Whole Genome and Transcriptome Amplification System



Enables whole genome and transcriptome sequencing from a single cell

Uses a single cell for the construction of a whole-genome and full-length mRNA transcriptome library



Provides industry-leading genomic coverage and resolution

Leverages a novel patented technology, primary template-directed amplification (PTA), to dramatically increase genomic capture and coverage to 97%^{1,2}



Superior transcriptome capture and coverage

Increases gene body coverage, representation across transcript sizes, and variant calling versus droplet-based RNA sequencing methods^{3,4}



Fits into established laboratory protocols

Compatible with various methods of single-cell singulation, sequencing platforms, and downstream applications, including whole exome and panel-based sequencing⁴



Scales to experiment size

Low-cost, automation-compatible, scalable approach with up to 384 reactions



Bioinformatics analysis included

Bioinformatics analysis and data visualization through **BaseJumper®** bioinformatics platform



References:

1. Gonzalez-Pena V, et al. Proc. Natl. Acad. Sci. U.S.A. 2021; 118 (24): e2024176118; doi: 10.1073/pnas.2024176118
2. Luquette L, et al. Nat Gen. 2022; 54: 1564–1571. doi: 10.1038/s41588-022-01180-2
3. Marks JR, et al. bioRxiv. 2023; doi: <https://doi.org/10.1101/2022.04.29.489440>
4. Data on file

Assay Performance

ResolveOME WGS DNA Performance	
Accuracy	99.99%
Sensitivity	96.65%
Specificity	99.99%
Allelic Balance	91.20%
Genomic Coverage	97.59%

ResolveOME WTS RNA Performance	
Protein Coding Genes	3451 ± 732
Concordance	0.97
Variance (CV)	32.9%

Figure 1: Assay Performance Characteristics. Analysis of NA12878 single cells prepared with ResolveOME Whole Genome Single-Cell Core Kit versus gold standard reference. WGS: whole genome sequencing, WTS: whole transcriptome sequencing.

A Revolution in Resolution From Each Cell

DNA

- Resolve SNV
- Resolve SV
- Resolve CNV
- Resolve Ploidy
- Resolve Exomes
- Resolve Genomes
- Resolve Panels
- Resolve Edits

DNA + RNA

- Resolve Transcriptomes
- Resolve Isoforms
- Resolve Cell ID
- Resolve Fusions

Resolve More.

ResolveServicesSM

Custom-built service projects, from singulating cells to figures. Services can include:

- Cell and/or nuclei sorting from fresh or frozen cells and tissues
- Whole genome amplification or whole genome and transcriptome amplification
- Exome capture or targeted protein analysis
- Library preparation for downstream applications
- Analysis using our bioinformatics platform, **BaseJumper[®]**, or bespoke computational analysis

Products

Codes	Product	Description
100956 100957	ResolveOME™ Whole Genome and Transcriptome Single-Cell Core Kit	PTA-based kit for whole genome and transcriptome amplification plus NGS library preparation from single cells. 96 or 384 reaction formats.
100954 100955	ResolveDNA® Whole Genome Single-Cell Core Kit	PTA-based kits for whole genome amplification plus NGS library preparation from single cells. 96 or 384 reaction formats.
100545	ResolveDNA® Whole Genome Amplification Kit	PTA-based kit for whole genome amplification from as little as 4 pg input DNA. Capable of 96 or 384 reactions.
100605	BaseJumper® Bioinformatics Platform	A platform for multiomic data analysis and visualization.
Early- Access	ResolveXOME™ Exome Capture Module	Exome capture module for use with ResolveDNA or ResolveOME Single-Cell Core Kits.

For a complete list of services, products, and pricing, email a member of our team, info@bioskryb.com

BioSkryb
GENOMICS

bioskryb.com

All data on file.
BIOSKRYB, RESOLVEDNA, RESOLVEOME, RESOLVEXOME, RESOLVESERVICES, and BASEJUMPER are trademarks of BioSkryb, Inc.
All other product names and trademarks are the property of their respective owners.
© 2024 BioSkryb, Inc. All Rights Reserved

